

Selectin Biosciences Inc.

Endo- β -Galactosidase

Recombinant from *Bacteroides fragilis*

Endo- β -Galactosidase (EC 3.2.1.103) cleaves internal β (1-4) galactose linkages in unbranched, repeating poly-N-acetylglucosamine structures. Sulfated structures such as keratan sulfate are also cleaved. Branching and/or fucosylation of the substrate may decrease or eliminate cleavage.

Endo- β -Galactosidase is useful for identifying and removing poly-N-acetylglucosamine structures on many biologically important glycoconjugates.

Product Code: GE 38

Specifications

Activity: 150 U/mg 15 U/mL

Storage: Store at 4°C. Do not freeze.

Formulation: The enzyme is provided as a sterile solution in 20 mM Tris-HCl pH 7.5.

Stability: Stable at least 12 months when stored properly. Several days exposure to ambient temperature will not reduce activity. Active for at least 5 days under reaction conditions.

Product Description

Molecular Weight

32,000 Daltons

Purity

Endo- β -Galactosidase is tested for contaminating protease as follows: 10 μ g of denatured BSA is incubated for 24 hours at 37°C

with 2 μ L of enzyme. SDS-PAGE analysis of the treated BSA shows no evidence of degradation.

The production strain of *E. coli* has been extensively tested and does not produce any detectable glycosidases.

Specificity: Internal β (1-4) galactose linkages in unbranched, repeating poly-N-acetylglucosamine [GlcNAc β (1-3)Gal β (1-4)]_n structures are the preferred substrate. Sulfated structures such as keratan sulfate are also cleaved. Branching and/or fucosylation of the substrate may decrease or eliminate cleavage. Sulfation of C-6 on galactose will block cleavage. Oligosaccharides of the neo-lacto group are cleaved at greatly reduced rates depending on the deviation from the preferred substrate. For example, Gal β (1-3)GlcNAc β (1-3)Gal β (1-4)Glc is cleaved at 5X10⁻⁵ the rate of keratan sulfate (see ref.4). Specificity is similar to the *Escherichia freundii* enzyme except that it is limited to cleaving N-acetylglucosamine extensions on tetraantennary structures of erythropoietin (see ref 5).

pH Range: Optimum: pH 5.8

Assay: One unit of Endo- β -Galactosidase is defined as the amount that will liberate one μ mole of reducing sugar per minute at 37°C and pH 5.8 from bovine corneal keratan sulfate.

Reagents

- 5X Reaction buffer is 250 mM sodium phosphate pH 5.8

Selectin Biosciences Inc.

Suggestions for Use:

Procedure for glycoproteins:

1. Add up to 100 µg of glycoprotein to a tube.
2. Add 4 ul 5X buffer and water to 19 µl.
3. Add 1 µl enzyme.
4. Incubate at 37°C for 2 hrs.

Procedure for oligosaccharides: Same as above except incubate from several hours to several days depending on the substrate. Add bovine serum albumen to 2 mg/ml to stabilize the protein during extended incubations.

References:

1. Scudder, P., Uemura, K., Doby, J., Fukuda, M.N. & Feizi, T. (1983) Isolation and characterization of an endo-β-galactosidase from *Bacteroides fragilis* **Biochem. J.** **213**, 485-494
2. Scudder, P., Hanfland, P., Uemura, K. & Feizi, T. (1984) Endo-β-galactosidases of *Bacteroides fragilis* and *Escherichia freundii* hydrolyze linear but not branched oligosaccharide domains of glycolipids of the neolacto series. **J. Biol. Chem.** **259**, 6586-6592.
3. Scudder, P., Tang, P.W., Hounsell, E.F., Lawson, A.M., Mehmet, H. & Feizi, T. (1986) Isolation and characterization of sulfated oligosaccharides released from bovine corneal keratan sulphate by the action of endo-β-galactosidase. **Eur. J. Biochem.** **157**, 365-373.
4. Murata, T., Hattori, T., Amarume, S., Koicki, A. & Usui, T. (2003) Kinetic

studies on endo-β-galactosidase by a novel colorimetric assay and synthesis of N-acetyllactosamine-repeating oligosaccharide β-glycosides using its transglycosylation activity. **Eur. J. Biochem** **270**, 3709-3719.

5. Hokke, C.H., Bergwerff, A.A., Van Dedem, D.W., Kamerling, J.P., and Vliegthart, J.F. (1995) Structural analysis of the N- and O-linked carbohydrate chains of recombinant human erythropoietin expressed in Chinese hamster ovary cells. Sialylation patterns and branch location of dimeric N-acetyllactosamine units. **Eur. J. Biochem.** **228**, 981-1008.

Selectin Biosciences Inc warrants that the above product conforms to the specifications described herein. Should the product fail for reasons other than through misuse Selectin Biosciences Inc. will, at its option, replace free of charge or refund the purchase price. This warranty is exclusive and Selectin Biosciences Inc. makes no other warrants, expressed or implied, including any implied conditions or warranties of merchantability or fitness for any particular purpose. Selectin Biosciences Inc. shall not be liable for any incidental, consequential or contingent damages.

This product is intended for in vitro research only.

REVISION 10/1/15

Selectin Biosciences Inc.