

# Selectin Biosciences Inc.

## $\alpha(2-3)$ Neuraminidase

$\alpha(2-3)$  Neuraminidase (N-acetylneuraminidase glycohydrolase EC 3.2.1.18) cleaves exclusively the non-reducing terminal  $\alpha(2-3)$  unbranched sialic acid residues from complex carbohydrates and glycoproteins. There is no detectable activity on  $\alpha(2-6)$  or  $\alpha(2-8)$  linkages or on branched  $\alpha(2-3)$  linkages (see Figure 1). To cleave all non-reducing terminal sialic acid residues including branched sialic acids (linked to an internal residue) from complex carbohydrates and glycoproteins, use  $\alpha(2-3,6,8,9)$  Neuraminidase.

$\alpha(2-3)$  Neuraminidase is isolated from a clone of *Streptococcus pneumonia* (formerly *Diplococcus pneumonia*). The enzyme has been extensively characterized using oligosaccharide standards.

$\alpha(2-3)$  Neuraminidase is useful for:

- Structural analysis of oligosaccharides
- Determining sialic acid linkage
- Glycoprotein desialylation
- Removing heterogeneity from glycoproteins

**Product Code:** GE 20

## Specifications

**Activity:**  $\geq 150$  U/mg,  $>5$  U/mL

**Storage:** Store at 4°C. Do not freeze.

**Formulation:** The enzyme is provided as a sterile-filtered solution in 50 mM Sodium Phosphate pH 7.5.

**Stability:** Stable at least 12 months when stored properly. Several days exposure to ambient temperatures will not reduce activity.

## Product Description

**Molecular weight:** ~75,000 Daltons

**Specificity:** Only non-reducing terminal unbranched  $\alpha(2-3)$  sialic acids (see Figure 1).

**Purity:** Each lot of  $\alpha(2-3,6)$  Neuraminidase is tested for contaminating protease as follows: 10  $\mu$ g of denatured BSA is incubated for 24 hours with 2  $\mu$ L of enzyme. SDS-PAGE analysis of the treated BSA shows no evidence of degradation.

The production host strain has been extensively tested and does not produce any detectable glycosidases.

**pH Range:** 50 mM sodium phosphate (pH 6.0) provides the optimal buffer for enzyme activity with 3'-sialyllactose, a standard substrate. If glycosidase treatment is performed at suboptimal pH because of glycoprotein solubility or activity requirements, expect some diminution in enzyme activity.

## Assay

One unit of  $\alpha(2-3)$  Neuraminidase is defined as the amount of enzyme required to produce 1  $\mu$ mole of methylumbelliferone in 1 minute at 37°C, pH 5.0 from MU-NANA [2'-(4-methylumbelliferyl)- $\alpha$ -D-N-acetylneuraminic acid].

## Reagents

- 5X reaction buffer 6.0 – 250 mM sodium phosphate, pH 6.0

# Selectin Biosciences Inc.

## Procedure for De-sialylation

1. Add up to 100 µg of glycoprotein or 1 nmol of oligosaccharide to tube.
2. Add water to a total volume of 14 µL
3. Add 4 µL 5X Reaction Buffer 5.0.
4. Add 2 µL of α(2-3) Neuraminidase.
5. Incubate at 37°C for 1 hour.

Desialylation may be monitored by SDS-PAGE if the size differential between native and desialylated protein is sufficient for detection.

## References

1. Corfield, A. P., H. Higa, J. C. Paulson and R. Schauer. The specificity of viral and bacterial sialidases for alpha(2-3) and alpha(2-6)-linked sialic acids in glycoproteins. **Biochim Biophys Acta** **744**:121-126 (1983).
2. Dwek, R. A., C. J. Edge, D. J. Harvey, M. R. Wormald and R. B. Parekh. Analysis of glycoprotein-associated oligosaccharides. **Ann Rev Biochem** **62**:65-100 (1993).
3. Kobata, A. Use of endo- and exoglycosidases for structural studies of glycoconjugates. **Anal Biochem** **100**:1-14 (1979).
4. Glasgow, L R., J. C. Paulson and R. L. Hill. Systematic purification of five glycosidases from Streptococcus pneumonia. **J. Biol Chem** **252**:8615-8623 (1977).

5. Kelly, R. T., D. Greiff and S. Farmer. Neuraminidase activity in Diplococcus pneumonia. **J Bacteriol** **91**:601-3 (1965).
6. Prime, S. J. Dearnley, A. M. Venton, R. B. Parekh and C. J. Edge. Oligosaccharide sequencing based on exo- and endoglycosidase digestion and liquid chromatographic analysis of the products. **J Chromatogr A** **720**:263-274 (1996).


Selectin Biosciences Inc warrants that the above product conforms to the specifications described herein. Should the product fail for reasons other than through misuse Selectin Biosciences Inc. will, at its option, replace free of charge or refund the purchase price. This warranty is exclusive and Selectin Biosciences Inc. makes no other warrants, expressed or implied, including any implied conditions or warranties of merchantability or fitness for any particular purpose. Selectin Biosciences Inc. shall not be liable for any incidental, consequential or contingent damages.

This product is intended for in vitro research only.

REVISION 5/22/12

# Selectin Biosciences Inc.

Figure 1 – Linkage specificities showing cleavable residues (in bold) for  $\alpha(2-3)$  Neuraminidase

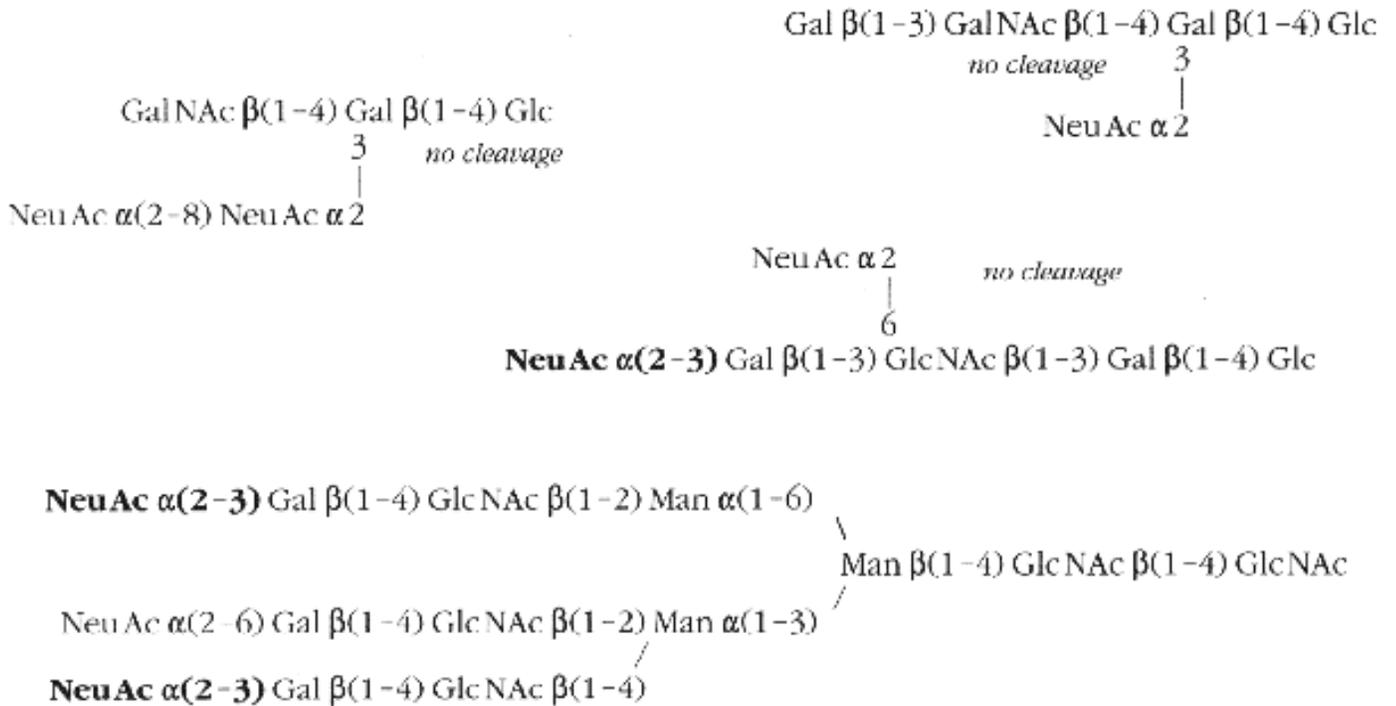

Gal = Galactose; Glc = Glucose; Man = Mannose; GalNAc = N-acetylgalactosamine; GlcNAc = N-acetylglucosamine; NeuAc = N-acetylneuraminic Acid (Sialic Acid)