

Selectin Biosciences Inc.

α (2-3, 6) Neuraminidase

α (2-3, 6) Neuraminidase (N-acetylneuraminidase, EC 3.2.1.18) cleaves all non-reducing terminal non-branched α (2-3)- and α (2-6) sialic acid residues from complex carbohydrates and glycoproteins. There is no detectable activity on α (2-8) or α (2-9) linkages or on branched α (2-3) or α (2-6) linkages. The relative cleavage rates for different linkages are: α (2-3) > α (2-6).

α (2-3, 6) Neuraminidase will not cleave branched sialic acids (linked to an internal residue). Use α (2-3, 6, 8, 9) Neuraminidase for α (2-8) or branched sialic acids. To cleave only non-reducing terminal α (2-3) unbranched sialic acid residues, use α (2-3) Neuraminidase.

α (2-3, 6) Neuraminidase is isolated from a clone of *Clostridium perfringens*. The enzyme has been extensively characterized using oligosaccharide standards.

α (2-3, 6) is useful for:

- Structural analysis of oligosaccharides
- Determining sialic acid linkage
- Glycoprotein deglycosylation
- Removing heterogeneity from glycoproteins

Product Code: GE 24

Specifications

Activity: ≥ 250 U/mg, ≥ 15 U/mL

Storage: Store at 4°C. Do not freeze.

Formulation: The enzyme is provided as a sterile-filtered solution in 20 mM Tris HCl, 25 mM NaCl pH 7.5.

Stability: Stable at least 12 months when stored properly. Several days exposure to ambient temperatures will not reduce activity.

Product Description

Molecular weight: ~41,000 Daltons

Specificity: Only non reducing terminal α (2-3) and α (2-6) unbranched sialic acids (Figure 1).

Purity: Each lot of α (2-3,6) Neuraminidase is tested for contaminating protease as follows: 10 μ g of denatured BSA is incubated for 24 hours with 2 μ L of enzyme. SDS-PAGE analysis of the treated BSA shows no evidence of degradation.

The production host strain has been extensively tested and does not produce any detectable glycosidases.

pH Range: 50 mM sodium phosphate (pH 6.0) provides the optimal buffer for enzyme activity with 3'-sialyllactose, a standard substrate. If glycosidase treatment is performed at suboptimal pH because of glycoprotein solubility or activity requirements, expect some diminution in enzyme activity.

Assay: One unit of α (2-3, 6) Neuraminidase is defined as the amount of enzyme required to produce 1 μ mole of methylumbelliferone in 1 minute at 37°C, pH 5 from MU-NANA [2'-(4-methylumbelliferyl)- α -D-N-acetylneuraminic acid].

Selectin Biosciences Inc.

Reagents

- 5X reaction buffer 6.0 – 250 mM sodium phosphate, pH 6.0

Suggestions for Use

Procedure for De-sialylation

Add up to 100 µg of glycoprotein or 1 nmol of oligosaccharide to tube

Add water to a total of 14 µL

Add 5 µL 5X Reaction Buffer 5.0

Add 2 µL of α(2-3,6) Neuraminidase

Incubate at 37°C for 1 hour.

Desialylation may be monitored by SDS-PAGE if the size differential between native and desialylated protein is sufficient for detection.

References

1. Corfield, A. P., H. Higa, J. C. Paulson and R. Schauer. The specificity of viral and bacterial sialidases for alpha(2-3) and alpha(2-6)-linked sialic acids in glycoproteins. **Biochim Biophys Acta** **744**:121-126 (1983).
2. Dwek, R. A. , C. J. Edge, D. J. Harvey, M. R. Wormald and R. B. Parekh. Analysis of glycoprotein-associated oligosaccharides. **Ann Rev Biochem** **62**:65-100 (1993).
3. Kobata, A. Use of endo- and exoglycosidases for structural studies of glycoconjugates. **Anal Biochem** **100**:1-14 (1979).
4. Prime, S. J. Dearnley, A. M. Venton, R. B. Parekh and C. J. Edge. Oligosaccharide sequencing based on exo- and endoglycosidase digestion and liquid chromatographic analysis of the products. **J Chromatogr A** **720**:263-274 (1996).
5. Roggentin, P, B. Rothe, F Lottspeich and R. Schauer. Cloning and sequencing of a Clostridium perfringens sialidase gene. **FEBS Lett** **238**:31-34 (Sep 1988).
6. Roggentin P., R. G. Kleineidam and R. Schauer. Diversity in the properties of two sialidase isoenzymes produced by Clostridium perfringens spp. **Biol chem Hoppe Seyler** **376**:569-575 (1995)


Selectin Biosciences Inc warrants that the above product conforms to the specifications described herein. Should the product fail for reasons other than through misuse Selectin Biosciences Inc. will, at its option, replace free of charge or refund the purchase price. This warranty is exclusive and Selectin Biosciences Inc. makes no other warrants, expressed or implied, including any implied conditions or warranties of merchantability or fitness for any particular purpose. Selectin Biosciences Inc. shall not be liable for any incidental, consequential or contingent damages.

This product is intended for in vitro research only.

REVISION 5/22/12

Selectin Biosciences Inc.

Figure 1 – Linkage specificities showing cleavable residues (in bold) for $\alpha(2-3, 6)$ Neuraminidase

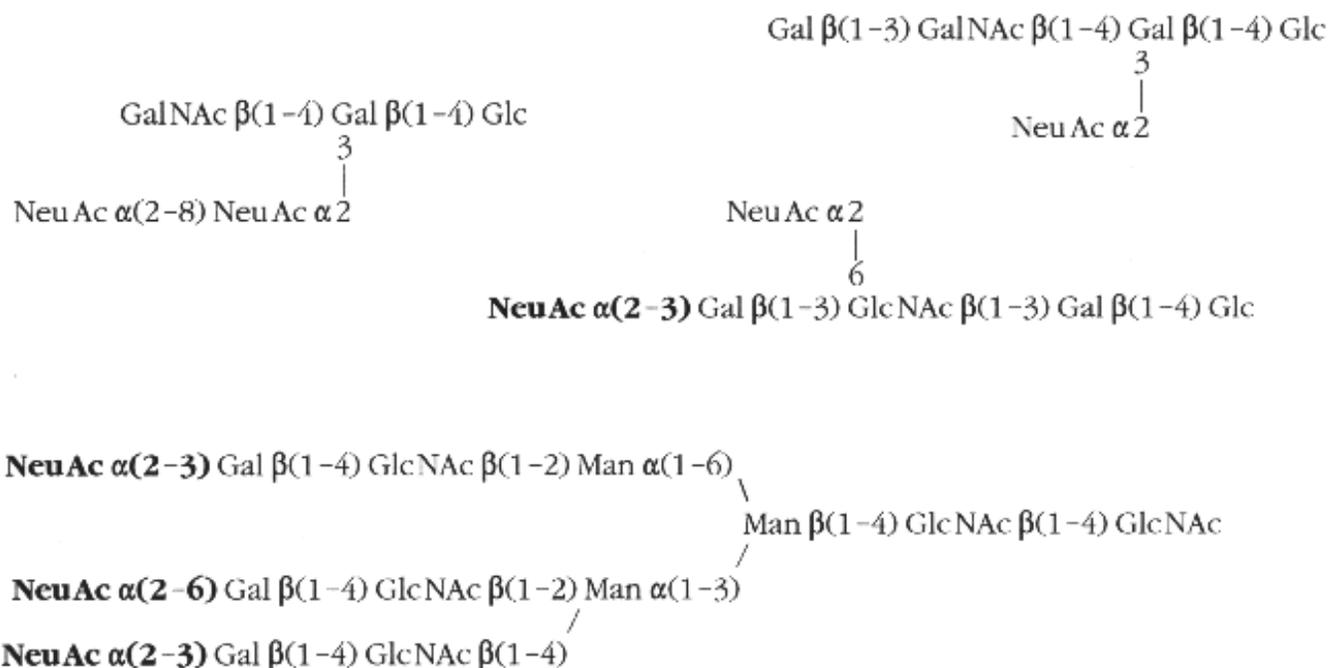

Gal = Galactose; Glc = Glucose; Man = Mannose; GalNAc = N-acetylgalactosamine; GlcNAc = N-acetylglucosamine; NeuAc = N-acetylneuraminic Acid (Sialic Acid)